DOUBLE VALVE 

DOUBLE VALVE GASKET 

CORA

CO.RA. ALTORNSCIO (LUCCA) ITALIA

DOUBLE VALVE

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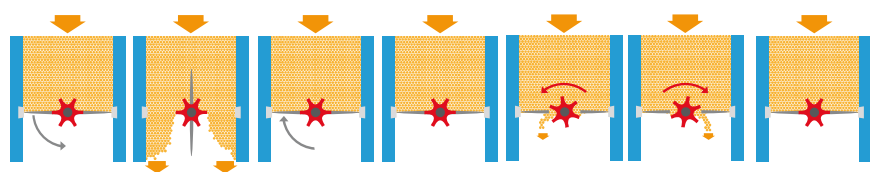
DOUBLE VALVE GASKET

The CO.RA. DOUBLE VALVE was conceived to satisfy control and dosing as well as packaging of loose powders during filling processes.

It consists of a butterfly with central rotor and is normally located on the bottom of machine process outlets such as driers or bins. This valve was designed to be inter-faced with an automatic and programmable weight system.

How the valve functions:

1. Heavy dosing:

The butterfly opens and discharges 95% of the pre-set gross weight. When the bulk weight limit is obtained the butterfly closes.

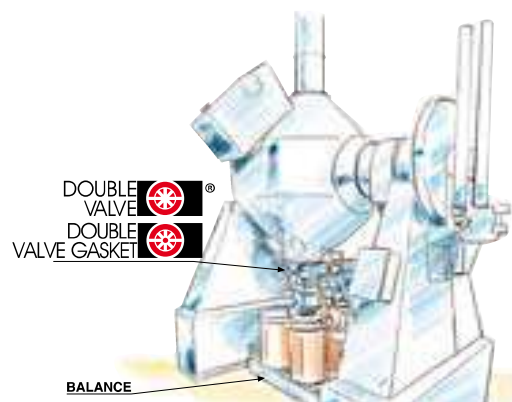
2. Fine dosing:

When the butterfly closes, the central rotor begins to trickle feed until the total weight of the product is obtained, with an accuracy which varies from ± 20 gr to ± 100 gr in compliance with characteristic of treated product. All is driven by 90° pneumatic actuator for the butterfly and 180° for the rotor.

The fine dosing pockets of the rotor can be customized in number and volume based on customer requirements. The most common application is the continuous liner dosing system, conceived by CO.RA., which allows the discharge of product inside a polyethylene liner as well as the filling of bags in a closed cycle, controlling the weight with great accuracy.








The DOUBLE VALVE GASKET is a variant of the Double Valve and has a seal between the butterfly and the dosing rotor. This solution allows the valve to work in applications with pressure up to 0.5 bar and with less product leakage during the discharge cycle.

Following numerous test we can declare that it is possible to use this type of valve in environments classified PBOEL 4.

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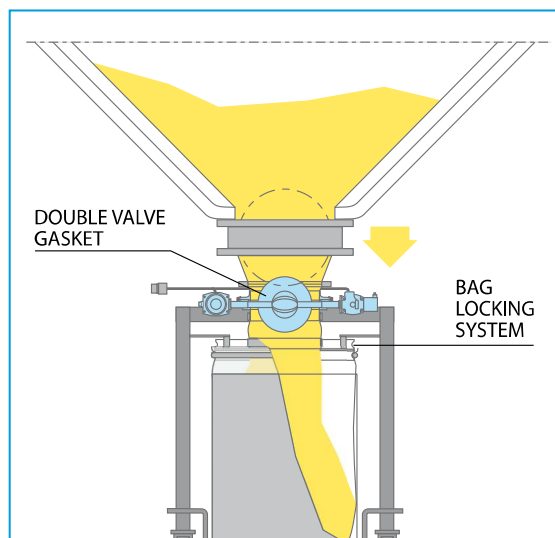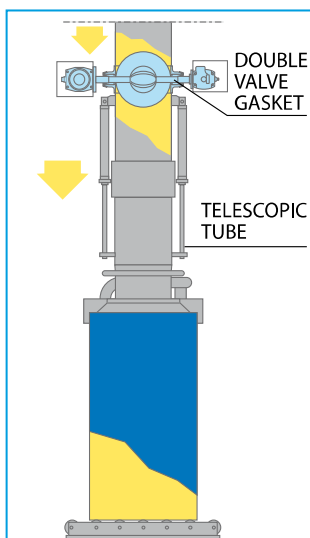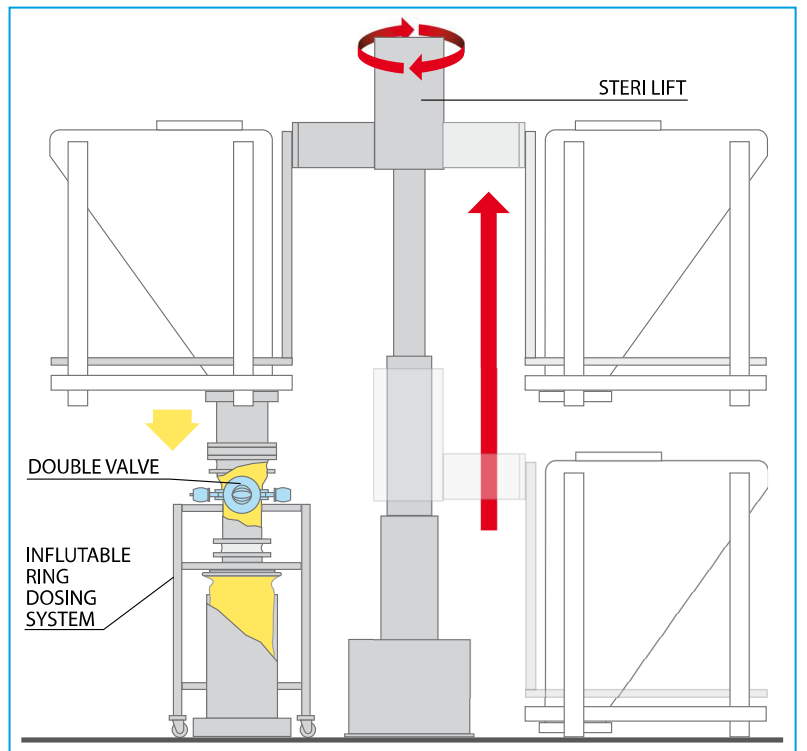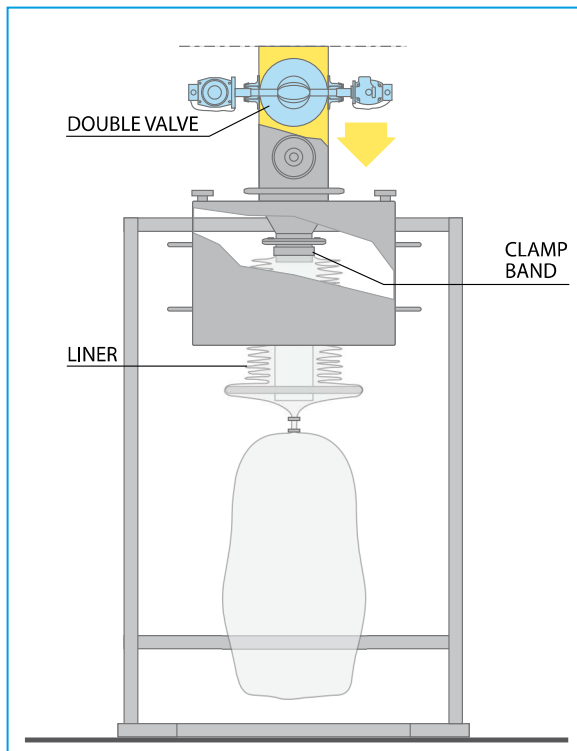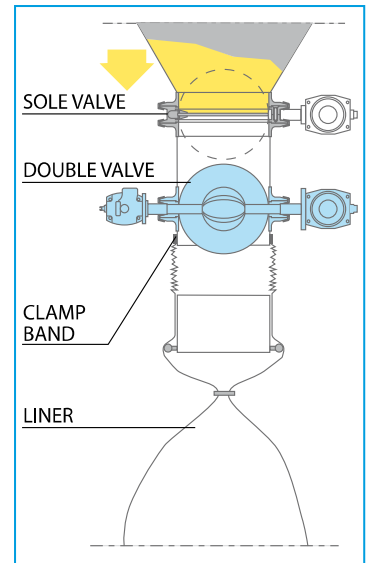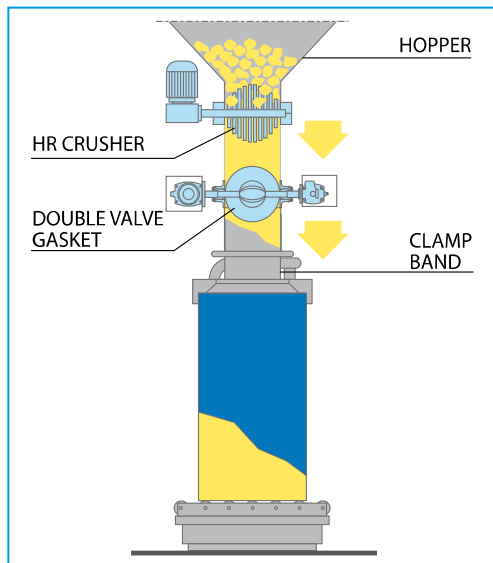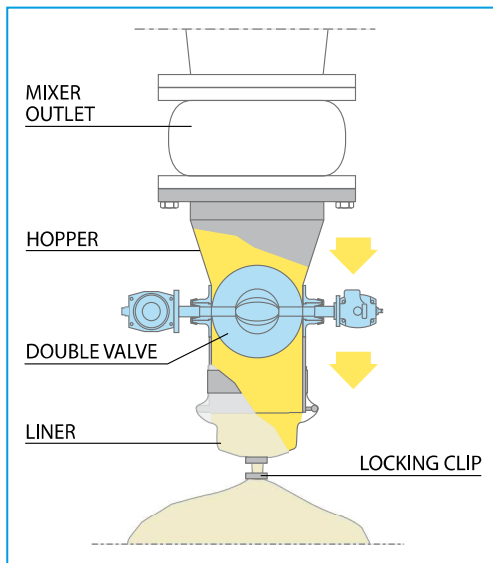For further information please refer to our "Pharmaceutical Process" brochure

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For further information please refer to our "Chemical Process" brochure

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- *Satin finish half-bodies in Aisi 316L (DIN 1.4404) mirror polished upon request, butterfly and rotor in mirror polished Aisi 316L (DIN 1.4404), other components in Aisi 304.*
- *Gaskets: platinum cured silicone, black silicone, VITON, EPDM, PTFE core with white or black silicone, silicone with stainless steel mesh*
- *Materials: stainless steel and silicone (certificates on request)*
- *Anodized aluminum pneumatic actuator, stainless steel on request*
- *Compressed Air: 6 bar*
- *NO O-ring on the shaft*
- *Easy to disassemble, inspect and clean*
- *Interchangeability with the SANITARY ROTARY FLOW (SRF)*
- *Very light weight and compact*
- *Ideal for granules and powders*
- *Manufactured according to G.M.P.*
- *Each valve bears the CO.RA. trademark and identification batch number (ex.: CO.RA. I/00/298/1)*

Advantages:

- *No lubricants*
- *Sterilizable in autoclave*
- *Minimum friction between butterfly and seal due to a reduced contact surface*

DIMENSIONI E AZIONAMENTI DOUBLE VALVE
DOUBLE VALVE DIMENSIONS AND ACTUATION

DIMENSIONI E AZIONAMENTI DOUBLE VALVE GASKET
DOUBLE VALVE GASKET DIMENSIONS AND ACTUATION

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On request it is possible to supply valves with WCWC - WCRC - RCRC half-bodies .

Attention: it is possible to supply a Ø 100 rotor for the Double Valve and Double Valve Gasket on request.

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The pneumatic actuator is the driving element that controls the movements of the butterfly and the rotor. The actuator for the control of the butterfly is normally double acting (D.E.) but a spring return (S.E.) actuator is available on request, while the actuator for the control of the rotor is always double acting (D.E.).

Air supply: filtered compressed air, better if lubricated, working pressure min. 5 bar/max. 6 bar.

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The Double Valve and Double Valve Gasket can be supplied with the following components on request:

- Actuator with stainless steel casing
- Namur 3/2, 5/2, 5/3 mono or bi-stable solenoid - standard or EX rated (Ex proof)
- Inductive proximity sensors, pneumatic and electric switches

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For more information regarding accessories please see page 6 of the Rotary Valve catalogue.

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THE DOUBLE VALVE AND DOUBLE VALVE GASKET IN THE DOSING SYSTEM

The DOSING SYSTEM allows the discharge and dosing of products such as powders and granules in bags, containers or drums that must be temporarily stored before shipping to the end-user, who could require high accuracy weight.

The DOUBLE VALVE and DOUBLE VALVE GASKET may be interfaced with the following Dosing Systems:

- Inflatable ring dosing system
- Continuous liner dosing system

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Rotore standard
Standard rotor

ROTOR AND DOSING ACCURACY

The dosing accuracy of the DOUBLE VALVE and DOUBLE VALVE GASKET is ± 50 grams with 90% of pharmaceutical products.

CO.RA. has found a solution to obtain higher dosing accuracy to meet customer requirements. The solution consists of customizing the volume and number of the rotor pockets so it can reach a dosing accuracy of ± 10 grams. In these cases it is necessary to "fine tune" the system by running tests using the customer's product. The demo tests can be carried out using either a DOUBLE VALVE or a DOUBLE VALVE GASKET that are available in our Technology Center.

The possibility to customize the number and volume of rotor pockets ensures the product a considerable performance flexibility, allowing the customer to optimize the process at any time simply by substituting the rotor.

Volume DOUBLE VALVE and DOUBLE VALVE GASKET rotor:

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*The DOUBLE VALVE and DOUBLE VALVE GASKET comply with Machinery Directive and are marked **CE**, they can also be supplied for ATEX zone. The valve markings can be as follows:*

II 1GD/2GD c 85°C (T6) 0°C Ta+40°
II 1D/2GD c 85°C (T6) 0°C Ta+40°
II 2GD/2GD c IIA 85°C (T6) 0°C Ta+40°
II 2GD/2GD c IIB 85°C (T6) 0°C Ta+40°

DOCUMENTATION AND CERTIFICATES

The documentation that can be supplied with the valve is:

- *3.1 material certificates for the parts in contact with the product and FDA and USP certificates for the gaskets*
- *Operation and Maintenance Manual*
- *Certificate of surface roughness*
- *Certificate of CE conformity*
- *FAT and SAT protocols*
- *IQ and OQ protocol*
- *Other documentation on request*

USEFUL INFORMATION

We suggest you to use the abbreviations indicated on page 8 (W, WC, ...) to easy fill in your P.O.

The components of the DOUBLE VALVE (DV) and DOUBLE VALVE GASKET (DVG) are interchangeable with an SRF of the same size.

Gasket available in several material and types are in accordance with FDA, for more information please ask for SOLE VALVE catalogues.

The DOUBLE VALVE and DOUBLE VALVE GASKET are designed and manufactured respecting customer requirements and to guarantee this, it is possible to run tests in our facilities with one of the DOUBLE VALVE or DOUBLE VALVE GASKET available in our Technology Center.

For further information please contact our Sales Dept. at +39 0583 20590 or e-mail: sales@coraitaly.net

REG. No. 5762/2

REG. No. IT-71044

Atex 2014/34/UE

CO.RA. S.r.l. - 55011 Spianate - Altopascio (Lucca) - Italy - Loc. Chiappini, 51
Tel. +39 0583 20590 r.a. - Fax +39 0583 20481 - info@coraitaly.net - www.coraitaly.net